


## An Introduction to Green Fluorescent Protein (GFP) | 1

---

Green fluorescent protein, more commonly known as GFP, was discovered in 1962. Since then, it has become one of the most widely exploited proteins for scientific research and has driven the development of many different microscopy-based applications. A limitation of GFP is that its fluorescence can be weakened or lost under certain fixation conditions. However, this problem can be resolved by using anti-GFP antibodies for signal enhancement.

## An Introduction to Green Fluorescent Protein (GFP) | 2

---

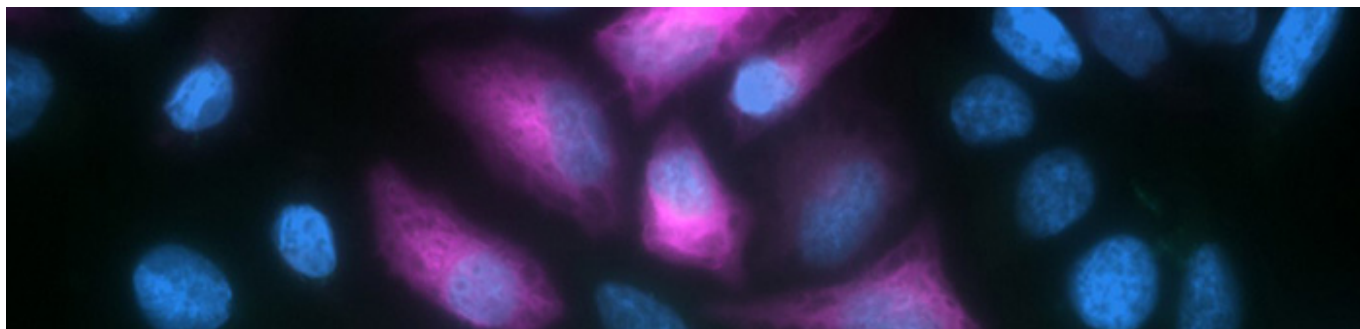

### A Brief History of GFP

## An Introduction to Green Fluorescent Protein (GFP) | 3

---

GFP was discovered by Osamu Shimomura in 1962 while investigating the natural bioluminescence of the jellyfish *Aequorea victoria*<sup>1</sup>. The aim of the original study was to identify the source of the blue light emitted by these organisms, which Shimomura found to be the Ca<sup>2+</sup> sensitive protein aequorin. However, during the purification of aequorin, a second protein that exhibited a bright green fluorescence was also isolated, albeit in only trace amounts.

In 1974, the green fluorescence of this second protein was shown to occur as a result of Förster resonance

## An Introduction to Green Fluorescent Protein (GFP) | 4

---

energy transfer (FRET) from aequorin<sup>2</sup>. But it was not until 1979 that the real breakthrough came, when Shimomura found the GFP chromophore to be part of the peptide chain, opening up the possibility of cloning the protein<sup>3</sup>. GFP was successfully cloned and sequenced by Prasher *et al.* in 1992<sup>4</sup>. Then, just two years later, Chalfie *et al.* expressed GFP in *Escherichia coli* and *Caenorhabditis elegans*<sup>5</sup>.

The structure of GFP was resolved by Tsien *et al.* in 1996, who showed the wild-type protein to consist of 238 amino acids (26.9kDa), folded to produce an  $\alpha$ -helix surrounded by a  $\beta$ -barrel<sup>6</sup>. The chromophore,

## An Introduction to Green Fluorescent Protein (GFP) | 5

---

located within the  $\alpha$ -helix (and thus protected by the  $\beta$ -barrel from fluorescent quenchers such as water, triplet oxygen, and photoisomerization), was found to be formed from serine 65, tyrosine 66, and glycine 67. GFP has major and minor excitation peaks at wavelengths of 395 nm and 475 nm, respectively, and emits fluorescence at 508nm, but only when the protein is correctly folded.

Tsien *et al.* subsequently engineered native GFP to be brighter and more photostable, as well as developed various GFP derivatives with different spectral characteristics<sup>7</sup>. Examples of these newer fluorescent

## An Introduction to Green Fluorescent Protein (GFP) | 6

---

proteins include enhanced GFP (eGFP), and cyan, blue, and yellow fluorescent proteins (CFP, BFP, and YFP, respectively). Shimomura, Chalfie, and Tsien were jointly awarded the Nobel Prize for Chemistry in 2008 for their discovery and development of the green fluorescent protein.

## An Introduction to Green Fluorescent Protein (GFP) | 7

---

GFP has been widely adopted for microscopy-based research as it enables the visualization of proteins of interest with no need for exogenous substrates or co-factors. The most common approach involves cloning the GFP gene downstream of a target gene in a cell line or model organism, whereby expression of the resultant fusion protein allows for monitoring target localization in either fixed or living samples.

A limitation of this method when working with fixed samples is that the fluorescence can be compromised by certain fixation conditions. One way of addressing this issue is to use anti-GFP antibodies to detect the

## An Introduction to Green Fluorescent Protein (GFP) | 8

---

expressed fusion protein. These can be either directly labeled with another fluorophore or paired with an appropriate secondary antibody for detection. An advantage of using secondary antibodies is that they can provide signal amplification.

For live-cell imaging applications, the time taken for GFP maturation prohibits the analysis of short-lived proteins, such as transcription factors, which are often degraded before the GFP produces a fluorescent signal. This problem is being overcome with the development of genetically encoded antibody-based



## An Introduction to Green Fluorescent Protein (GFP) | 9

---

probes such as antigen-binding fragments (Fabs), single-chain variable fragments (scFvs), and camelid nanobodies that are pre-labeled with mature GFP or other fluorophores<sup>8,9</sup>.

Another important application for GFP is its use for co-localization studies. Here, one protein is labeled with GFP while another is labeled with a different fluorophore such that the emission spectrum of one of the fluorophores overlaps with the excitation spectrum of the other. By illuminating the sample with a laser that excites just the first fluorophore, a signal will only be produced if the two proteins are in close enough

proximity for FRET to take place.

GFP has also been exploited for studies based on fluorescence recovery after photobleaching (FRAP) and fluorescence loss in photobleaching (FLIP). These techniques involve photobleaching a region of interest within a cell, often an area of the cell membrane, before monitoring the diffusion of unbleached GFP back into the area in question (FRAP) or tracking the decrease in fluorescence outside of the region of interest as the bleached proteins disseminate (FLIP).

## Supporting your research

GFP and its derivatives represent indispensable tools for scientific research that underpin both established and emerging applications. Jackson ImmunoResearch has recently launched a new anti-GFP antibody, which is available conjugated to a range of conjugates, including Alexa Fluor® dyes, or they can be paired with our high-quality anti-rabbit secondaries to amplify GFP detection.

[Learn more about Anti-GFP Antibodies](#)

**References:**

- [SHIMOMURA O, JOHNSON FH, SAIGA Y. Extraction, purification and properties of aequorin, a bioluminescent protein from the luminous hydromedusan, Aequorea. \*J Cell Comp Physiol.\* 1962;59:223-239. doi:10.1002/jcp.1030590302 <https://pubmed.ncbi.nlm.nih.gov/13911999/>](#)
- [Shimomura O, Johnson FH, Morise H. Mechanism of the luminescent intramolecular reaction of aequorin. \*Biochemistry.\* 1974;13\(16\):3278-3286. doi:10.1021/bi00713a016 <https://pubmed.ncbi.nlm.nih.gov/4152180/>](#)
- [O. Shimomura, Structure of the chromophore of Aequorea green fluorescent protein, \*FEBS Letters\*, Volume 104, Issue 2, 1979, Pages 220-222, <https://www.sciencedirect.com/science/article/pii/0014579379808182>](#)
- [Prasher DC, Eckenrode VK, Ward WW, Prendergast FG, Cormier MJ. Primary structure of the Aequorea victoria green-fluorescent protein. \*Gene.\* 1992;111\(2\):229-233. doi:10.1016/0378-1119\(92\)90691-h <https://pubmed.ncbi.nlm.nih.gov/1347277/>](#)
- [Chalfie M, Tu Y, Euskirchen G, Ward WW, Prasher DC. Green fluorescent protein as a marker for gene expression. \*Science.\* 1994;263\(5148\):802-805. doi:10.1126/science.8303295 <https://pubmed.ncbi.nlm.nih.gov/8303295/>](#)
- [Ormö M, Cubitt AB, Kallio K, Gross LA, Tsien RY, Remington SJ. Crystal structure of the Aequorea victoria green fluorescent protein. \*Science.\* 1996;273\(5280\):1392-1395. doi:10.1126/science.273.5280.1392 <https://pubmed.ncbi.nlm.nih.gov/8703075/>](#)
- [Tsien RY. The green fluorescent protein. \*Annu Rev Biochem.\* 1998;67:509-544. doi:10.1146/annurev.biochem.67.1.509 <https://pubmed.ncbi.nlm.nih.gov/9759496/>](#)
- [Lyon K, Stasevich TJ. Imaging Translational and Post-Translational Gene Regulatory Dynamics in Living Cells with Antibody-Based Probes. \*Trends Genet.\* 2017;33\(5\):322-335. doi:10.1016/j.tig.2017.02.003 <https://pubmed.ncbi.nlm.nih.gov/28359585/>](#)
- [Zhao N, Kamijo K, Fox PD, et al. A genetically encoded probe for imaging nascent and mature HA-tagged proteins in vivo. \*Nat Commun.\* 2019;10\(1\):2947. Published 2019 Jul 3. doi:10.1038/s41467-019-10846-1 <https://pubmed.ncbi.nlm.nih.gov/31270320/>](#)

**Learn more:**[Indirect and direct Western blotting](#)[Chemiluminescence western blotting](#)[An Introduction to Expansion Microscopy](#)**Do more:**[Exhibition Schedule](#)[Western blotting guide](#)[ELISA guide](#)